

# THz-Raman – Accessing molecular structure with Raman spectroscopy for enhanced chemical identification, analysis and monitoring

Randy A. Heyler<sup>a</sup>, James T.A. Carriere<sup>a</sup>, and Frank Havermeyer<sup>a</sup>

<sup>a</sup>Ondax Inc. 850 E. Duarte Road, Monrovia, CA 91016

## ABSTRACT

Structural analysis via spectroscopic measurement of rotational and vibrational modes is of increasing interest for many applications, since these spectra can reveal unique and important structural and behavioral information about a wide range of materials. However these modes correspond to very low frequency ( $\sim 5\text{cm}^{-1}$  -  $200\text{cm}^{-1}$ , or 150 GHz-6 THz) emissions, which have been traditionally difficult and/or expensive to access through conventional Raman and Terahertz spectroscopy techniques. We report on a new, inexpensive, and highly efficient approach to gathering ultra-low-frequency Stokes and anti-Stokes Raman spectra (referred to as “THz-Raman”) on a broad range of materials, opening potential new applications and analytical tools for chemical and trace detection, identification, and forensics analysis. Results are presented on explosives, pharmaceuticals, and common elements that show strong THz-Raman spectra, leading to clear discrimination of polymorphs, and improved sensitivity and reliability for chemical identification.

**Keywords:** Raman, Low-frequency Raman, THz spectroscopy, Polymorph Identification, Explosives Detection, Chemical Identification, Forensics, Volume Holographic Notch Filters

## 1. INTRODUCTION

Spectroscopic analysis of the molecular and intermolecular structure of a substance is of increasing importance for many applications, since these spectra can reveal unique information about the physical properties and behavior of an element or compound. Consider carbon, which can take the form of graphite, diamond, graphene, “Buckyball” or nanotube, among others, all of which behave quite differently. (Such substances are known generally as polymorphs, or in the case of a single element, an allotrope). Identification of polymorphs is critical in many applications, including formulation and analysis of pharmaceuticals, chemicals and polymers. And structural information can improve the sensitivity and accuracy of explosives detection, while revealing important clues about preparatory methods, manufacture techniques, or sources of origin during forensic analysis. However, getting access to molecular structural information via spectroscopy has typically been either difficult, expensive, and/or tethered to the laboratory.


Raman spectroscopy is a “scattering” spectroscopic technique that has been widely adopted as an economical and effective means of optically sensing and analyzing the chemical composition (or “fingerprint”) of a substance. The corresponding “fingerprint region” of the Raman spectrum ranges from about  $200\text{cm}^{-1}$  to  $2,000\text{cm}^{-1}$ . However, it has long been known that “low-frequency” Raman (defined as a Raman shift of  $\sim 5\text{cm}^{-1}$  to  $200\text{cm}^{-1}$ , or 150GHz-6THz in the electromagnetic spectrum) can yield rich information regarding the vibrational energies associated with molecular and intermolecular structures and features<sup>1</sup>. Unfortunately, most Raman systems – though widely adopted in a broad range of laboratory and field uses – remain limited to capturing only chemical fingerprint information, primarily due to the use of edge or notch filters that also remove most, if not all of these low frequency signals, including the entire anti-Stokes region. Some higher-end Raman systems can be configured to access the low-frequency regime, but these are mostly large and expensive laboratory instruments, and often have very limited throughput. Thus, most Raman systems leave out a vast region of important spectral information that could be used to improve sensitivity and accuracy, as well as yield insights into the structure of the materials they measure.

Alternatively, terahertz (THz) spectroscopy is an “absorption” technique that can be used to measure these structural modes by directly emitting radiation in the 300GHz to 6THz frequency range ( $10\text{cm}^{-1}$  to  $200\text{cm}^{-1}$  region), and detecting the corresponding absorption spectrum. THz systems have the additional benefit of being able to more deeply penetrate

a material or “see through” external layers (e.g. clothes, packaging, etc.) to capture the signal. Unfortunately, these systems rely on expensive laser sources, and limits on detector performance, availability and expense have constrained the potential sensitivity, resolution, and/or economics of using this technique. Plus, their fairly narrow spectral range (only up to 3 to 6 THz) limits their ability to do complete and reliable chemical identification for many materials.

“THz-Raman” refers to a new technique that extends the range of Raman spectroscopy from the fingerprint region into the Terahertz regime, adding important insight to molecular and intermolecular structure to the data about chemical composition (Figure 1). It has been shown that low-frequency Raman/THz-regime spectra can be used to greatly improve differentiation and analysis of both the structure and chemistry of materials, thereby improving accuracy, sensitivity, and scientific or forensic analysis<sup>2</sup>. Detection, identification, and forensic analysis of many substances, including explosives, drugs, pharmaceuticals, biological tissues, polymers, and hazardous materials, can all benefit from this extended spectral information.

We report here on a new, compact, efficient and affordable way to *simultaneously* deliver both chemical and structural analysis to a wide range of applications. An integrated, compact THz-Raman system is demonstrated utilizing a single stage spectrometer to show both Stokes and anti-Stokes measurements down to  $<10\text{cm}^{-1}$  on a wide variety of materials, including but not limited to explosives, pharmaceuticals, hazardous chemicals and biological specimens. These all are shown to exhibit characteristically different low frequency/THz peaks that can be exploited for chemical analysis, identification, monitoring, forensics, and contamination applications.

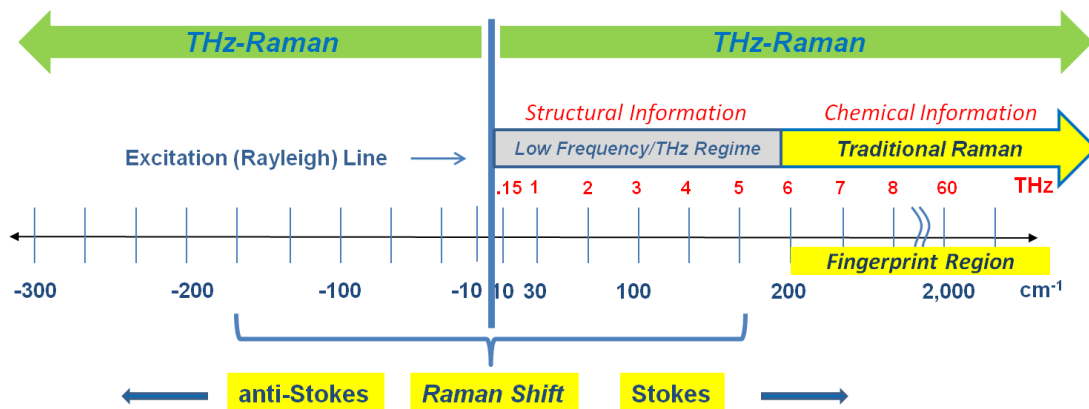


**Figure 1.** THz-Raman combines the entire low-frequency/THz regime, the traditional fingerprint region, and anti-Stokes spectral ranges, yielding additional structural information and improving chemical detection sensitivity.

## 2. THZ-RAMAN SPECTROSCOPY


As mentioned above, “THz-Raman” refers to the simultaneous capture of both Stokes and anti-Stokes shifts in the ultra-low frequency region (from  $\pm 5\text{cm}^{-1}$  to  $200\text{cm}^{-1}$ ). (THz-Raman systems also capture the “fingerprint” region from  $200\text{cm}^{-1}$  to  $>2,000\text{cm}^{-1}$ ). The low-frequency regions (for both Stokes and anti-Stokes shifts) are particularly difficult to resolve, as they reside extremely close to the Rayleigh excitation wavelength, which needs to be attenuated by at least OD 8 in order to see any Raman signals. (Raman scattering is a very weak phenomenon, with only about 1 in every  $10^8$  photons exhibiting a Raman shift). Most traditional Raman systems currently utilize thin-film edge filters, which end up completely removing both the Rayleigh excitation *and* the entire anti-Stokes region, cutting off all signals below about  $200\text{cm}^{-1}$  from the Rayleigh line. And most commercial notch filters, which will allow some of the anti-Stokes signal to come through, will still block all signals below about  $\pm 150\text{cm}^{-1}$ . Getting *both* high Rayleigh attenuation *and* resolution close to the excitation line has only been possible by using a multi-stage spectrometer system, which (1) is large, bulky, and expensive, (2) requires very precise alignment and set-up and (3) significantly reduces the overall throughput of the already weak Raman signal.

Recent advances in volume holographic grating (VHG) filter technology<sup>3,4,5</sup> have enabled the manufacture of exceptionally narrow bandwidth notch filters with very high throughput (Figure 2). Each VHG filter has a notch profile that is designed to diffract only one specific wavelength matching the laser and transmit all other wavelengths. The ultra-narrow transition bandwidth of these filters enables extremely high attenuation of the laser wavelength (>OD 4 for each filter), while maintaining very high transmission of nearby Raman signals beyond  $\sim 5\text{cm}^{-1}$  (Figure 2).

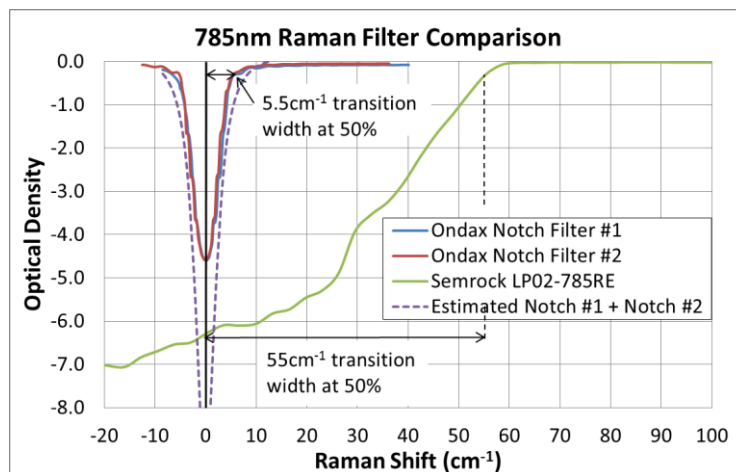

**Figure 2.** A VHG notch filter has an approximately 10x narrower transition width than a thin-film edge filter, enabling very high Rayleigh attenuation without blocking low-frequency or anti-Stokes Raman signals.

This has led to systems that are capable of rapid acquisition of high-quality, ultra-low frequency Raman spectra in the 5-200  $\text{cm}^{-1}$  region. These systems are based on a stable wavelength laser source, a series of VHG filters and a single stage spectrograph (Figure 3). This combination of strong Rayleigh attenuation and high broadband transmission enable the system to simultaneously capture both the intense low-frequency Stokes and anti-Stokes Raman bands and the “fingerprint region” transitions, greatly simplifying the overall system and reducing size and cost, while improving the sensitivity and reliability of using Raman for chemical identification and other applications.

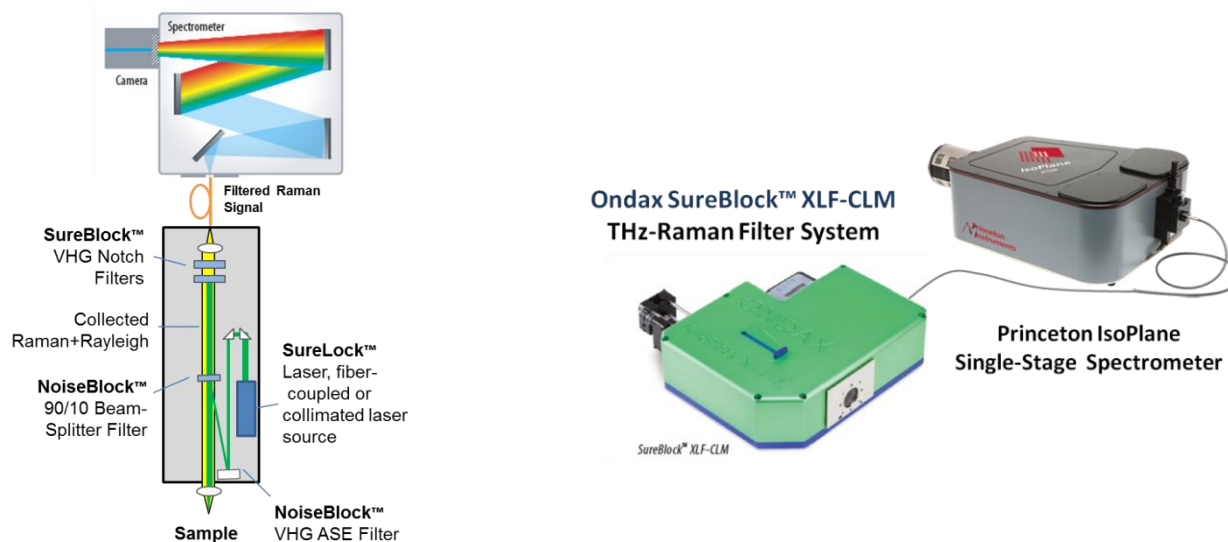

**Figure 3.** A THz-Raman system is comprised of a compact, confocal optical system that includes single-frequency laser source, ASE filters, a high-throughput beamsplitter, sample collection optics, and multiple cascaded VHG notch filters. The filtered Raman signal is fed into a single-stage spectrometer for readout. (A schematic of the system is depicted on the left, and a photograph of system is to the right).

### 3. THZ-RAMAN SYSTEM ARCHITECTURE

Experimental measurements were taken with the confocal THz-Raman system shown in Figure 3. The system was comprised of a single mode 785nm wavelength stabilized diode laser (SureLock™ LM series, Ondax, Inc.) and a series of ultra-narrowband VHG filters that were spectrally matched to the laser output wavelength. Two VHG ASE suppression filters (NoiseBlock™, Ondax, Inc.) were used to remove amplified spontaneous emission (ASE) from the laser. (ASE is often on the same order of magnitude or larger than the Raman signals, which can reduce SNR or swamp low frequency signals if it is not attenuated). A dichroic 90/10 VHG beamsplitter filter (NoiseBlock™, Ondax, Inc.) redirects the laser towards the sample, where a 10X objective lens focuses the laser onto the sample and collects the back-scattered light. The 90/10 beamsplitter then reflects 90% of the Rayleigh scatter back towards the laser while transmitting all of the Raman shifted signals. (The dichroic nature of the 90/10 beamsplitter results in an almost 4X improvement in collected Raman signal compared to a broadband 50/50 beamsplitter). Two ultra-narrowband VHG notch filters (SureBlock™, Ondax, Inc.), each having optical density >4.0, then further attenuate the collected Rayleigh scattered light while transmitting the Raman signals with an estimated system transmission efficiency of >80%. The filtered signal was focused into a 25µm core diameter, 0.1NA step index fiber (HPSC25, ThorLabs) and connected to a high-resolution, high-throughput single stage spectrometer (IsoPlane series, Princeton Instruments) 0.3m imaging spectrometer. It was equipped with a 1200 lines/mm grating and a 1340x400 imaging array (Pixis model 400BR with eXcelon, Princeton Instruments) with 20 x 20 µm pixel size and 98% peak quantum efficiency to ensure maximum signal collection and ~1.25cm<sup>-1</sup> resolution; appropriately matched for analysis of the 5-200cm<sup>-1</sup> frequency range. (Note: the range, resolution, and noise level of the overall system is ultimately determined by the specifications of the spectrometer, including grating, detector, and size). As indicated in the results section, some measurements were taken with a similar system configured for 532nm operation, using a diode-pumped solid state (DPSS) 532nm laser (GEM, Laser Quantum) and matching 532nm ASE, beamsplitter and notch filters. (These systems can be configured for 488, 514, 532, 633, 785 or 830nm operation).

The entire laser and filter assembly is extremely compact – approximately the size of a ream of notebook paper – and the fiber coupling allows for remote location of the spectrometer (we have demonstrated at least 15m separation without signal loss or interference from fluorescence). The low power requirements of the laser also enable the filter system to be battery operated if needed. Finally, it is important to note that the ultra-narrow bandwidth of the VHG filters (<0.1nm) require the use of either a wavelength-stabilized diode laser, narrow-band DPSS laser, or gas laser. Normal diode lasers tend to mode hop, shifting the laser wavelength outside the blocking range of the filters and resulting in either reduced Rayleigh suppression or a complete loss of signal in the described configuration.


### 4. EXPERIMENTAL RESULTS

#### 4.1 Polymorph Identification


As noted above, one key application of chemical structural analysis is the identification of polymorphs (defined as solid materials that exist in more than one form or crystal structure). For example, in the pharmaceutical industry many compounds exhibit polymorphism, and each form may exhibit significantly different therapeutic efficacy and bio-availability. Therefore only one form of the compound is normally approved for medical use. In the petrochemical industry, the form of the polymer chain can strongly affect the physical properties of the material. Even some explosives exhibit polymorphism. So it is often critical that the polymorphic form can be clearly and reliably identified during formulation, quality assurance, or analysis.

One good example is Carbamazepine (CBZ), an anticonvulsant and mood-stabilizing drug commonly prescribed for treatment of epilepsy and bipolar disorder. It has four different polymorphic forms that have been well characterized in the literature<sup>6,7,8,9,10</sup> with form 3 being the active pharmaceutical ingredient (API) approved for medical use. We obtained pure samples of form 2 and form 3 as well as the hydrated form, and measured the entire Raman spectra (Figures 4a and 4b). As expected, since the molecules have the same chemical composition, the fingerprint region signals are quite similar. However, the different structural forms of the polymorphs are clearly evident in the low frequency/THz regime, and are actually much stronger than those in the fingerprint region. Figure 4b shows expanded details of the various polymorphs and Hydrate in this region. Note that the anti-Stokes signals (which are symmetrical to

the Stokes signals but lower in intensity) can also be used to clearly validate the low-frequency measurements, as well as providing additional information that can be used to boost the detection capabilities of the system.

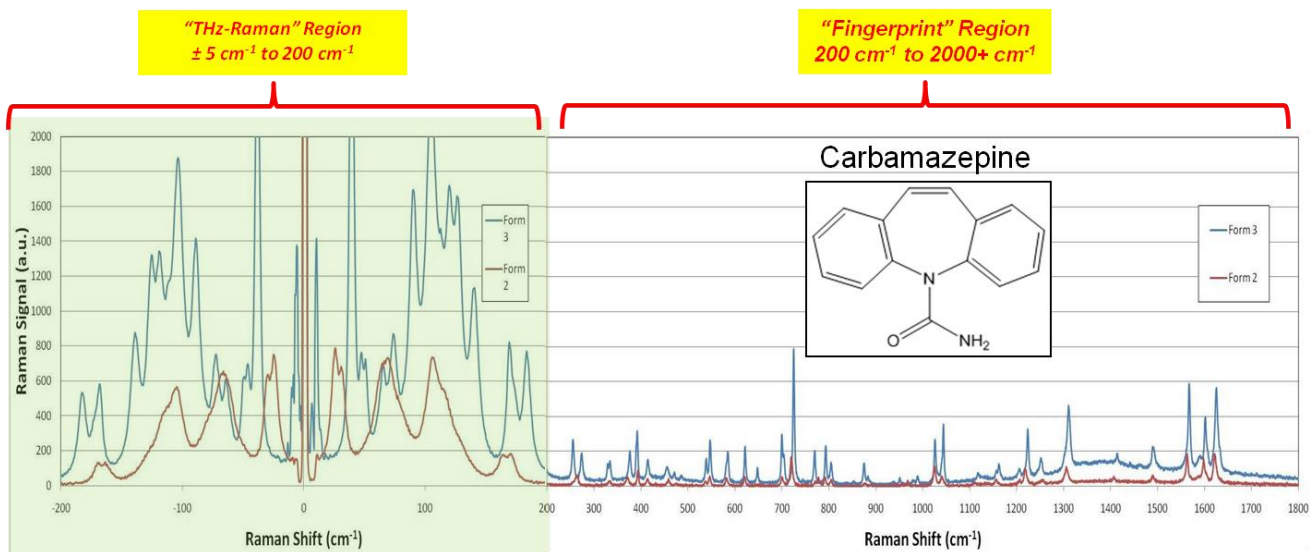


**Figure 4a.** The THz-Raman spectra of two polymorphs of Carbamazepine (CBZ), showing extremely strong and differentiated low-frequency signals when compared to the fingerprint region. (Data taken at 785nm).

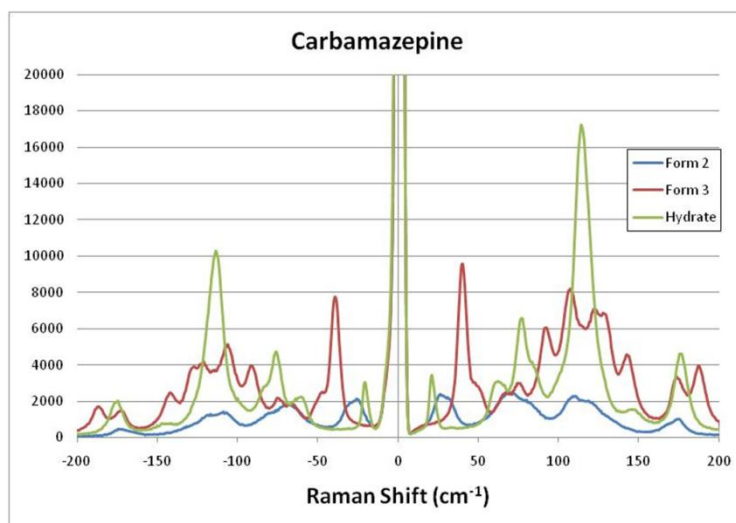

**Figure 4b.** A close-up of the low-frequency Stokes and anti-Stokes spectra of Form 2, Form 3 and the Hydrate of Carbamazepine (CBZ). (Data taken at 785nm).

When single elements such as Carbon exhibit polymorphism, they are called allotropes. An excellent example is Sulfur, which forms over 30 different allotropes<sup>11</sup>. The most common and easiest to produce are forms  $\alpha$ ,  $\beta$  and  $\lambda$ . We placed a sample of  $\alpha$  sulfur on a microscope slide and heated it while measuring the Raman spectra. When the sample temperature was increased above 95.2°C, the form changed from  $\alpha$  to  $\beta$ . Further increasing the temperature above the melting point at 115.2°C resulted in a second form change to  $\lambda$ . Figure 5 plots the corresponding Raman spectra for each form. While there is a change in magnitude of the peaks in the Raman fingerprint region for each form, there is no obvious shift in the position of the peaks. By comparison, the ultra-low-frequency peaks change dramatically from one form to another in both magnitude and location, enabling clear differentiation of the allotropes.

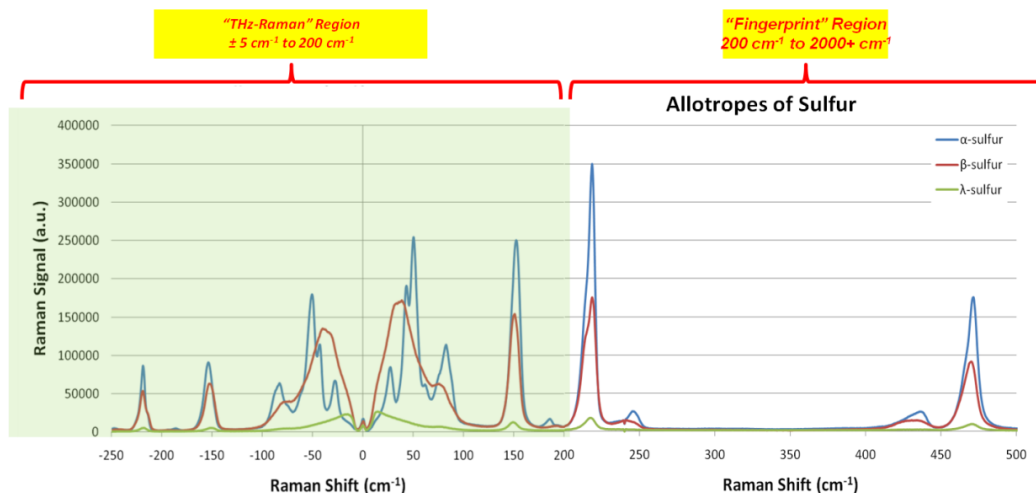


**Figure 5.** Allotropes of Sulfur. Sulfur transitions from  $\alpha$  to  $\beta$  form at 95.2 °C, and then to  $\lambda$  at 115.2 °C. (Data taken at 785nm).

## 4.2 Explosives and Drug Detection

Raman has been increasingly used for identification of explosives and illicit drugs in the security and defense field<sup>12</sup>. Portable, hand-held systems have now made this a ubiquitous tool for first-responders and military personnel around the world. However, there still remain some limitations to sensitivity and/or occasional false positives. THz-Raman offers the promise of not only improving overall sensitivity, but also providing a “forensic signature” that can reveal details about the structure, preparatory methods and other ingredients/contaminants that could help identify a source of origin of the material.

Ammonium Nitrate ( $\text{NH}_4\text{NO}_3$ ) is a common material used in homemade explosives (HME), specifically the Oklahoma City bombing in 1995 and also cited as the source material in the West, Texas blast in April, 2013. Figure 6 shows the Raman spectra of Ammonium Nitrate. Once again, notice the very strong, broad signals in the low frequency/THz-Raman region. Importantly, these signals will be even more pronounced (compared to the sharp spectra at higher frequencies) when observed with a lower-resolution system, since the signal is proportional to the total integrated area under the spectrum. Sharp, high peaks represent only a relatively small amount of photons, and while easily detectable in a high-resolution scan, they can nearly disappear when measured with a lower-resolution system, such as those commonly used in handheld systems. The broad peaks in the low frequencies will remain apparent under all conditions.

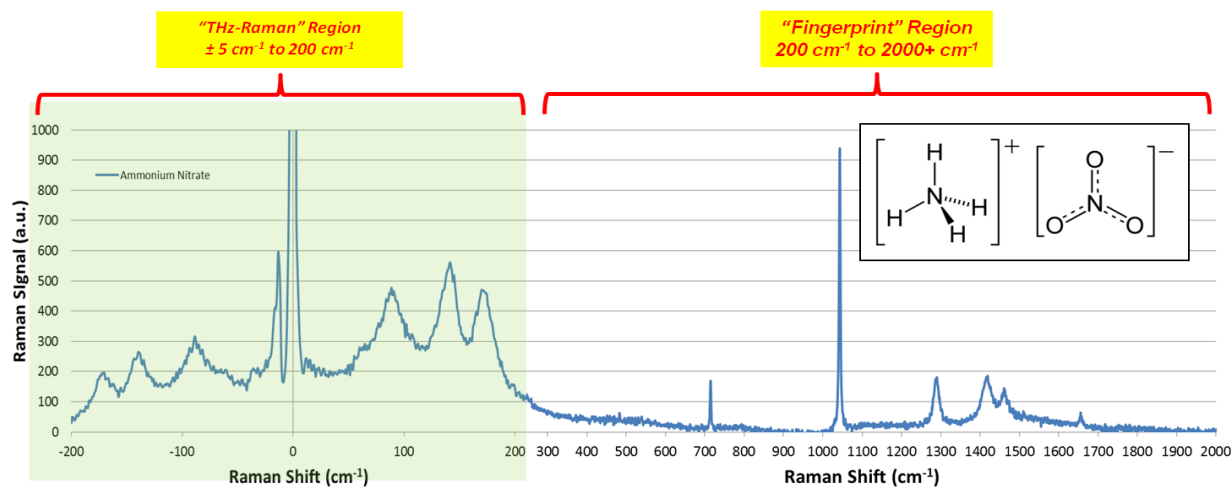

**Figure 6.** Ammonium Nitrate. The strong, broad peaks in the low-frequency/THz-Raman significantly boost overall signal and will remain observable even with a low-resolution scan. (Data taken at 532nm).

## 5. CONCLUSIONS

The addition of low-frequency/THz-Raman signals to traditional Raman spectra can boost the sensitivity and significantly enhance the forensic value of Raman spectroscopy for a wide range of applications, including polymorph identification, explosives trace detection, polymer and industrial chemical development and manufacture, material science, and forensics analysis. VHG-enabled THz-Raman systems are compact, efficient, and add a wealth of information regarding the physical structure of the molecule in addition to chemical composition. These new platforms greatly simplify the use and lower the cost of exploiting the many benefits of low-frequency Raman spectroscopy.

## 6. ACKNOWLEDGEMENTS

We would like to thank Princeton Instruments for providing the spectrometer and their ongoing support of our application development work.

---

<sup>1</sup> Hédoux, A et al “Using the low-frequency Raman spectroscopy to analyze the crystallization of amorphous Indomethacin,” European Journal of Pharmaceutical Sciences 38 156–164 (2009)

<sup>2</sup> Hédoux, A. et. al “The contribution of Raman spectroscopy to the analysis of phase transformations in pharmaceutical compounds” International Journal of Pharmaceutics 417 17– 31(2011)

<sup>3</sup> US Patents 7,986,407 and 8,184,285

<sup>4</sup> Moser, C., Havermeier, F., “Ultra-narrow band tunable laserline notch filter,” Applied Physics B: Lasers and Optics 95(3), pp. 597-601 (2009).

<sup>5</sup> Moser, et al “Compact Raman spectrometer system for low frequency spectroscopy,” 2010, Proc. SPIE, Vol.7598 (2010)

<sup>6</sup> Pohlmann H, Gulde C, Jahn R, Pfeifer S.”Polymorphism, particle-size and blood level values of carbamazepine,” Pharmazie 30:709–711 (1975)

<sup>7</sup> Lefebvre C, et al. “Polymorphic transitions of carbamazepine during grinding and compression,.) Drug Dev Ind Pharm 12:1913–1927.

<sup>8</sup> Rustichelli C, et al “Solid-state study of polymorphic drugs: Carbamazepine,.) J Pharm Biomed Anal 23:41–54.

<sup>9</sup> Lang M. et al “Form IV of carbamazepine” J Pharm Sci 91:1186–1190.

<sup>10</sup> Day et. al. “Understanding the Influence of Polymorphism on Phonon Spectra: Lattice Dynamics Calculations and Terahertz Spectroscopy of Carbamazepine,” J. Phys. Chem. B, 110, 447-456 (2006)

<sup>11</sup> Steudel, Ralf; Eckert, Bodo "Solid Sulfur Allotropes Sulfur Allotropes," Topics in Current Chemistry 230: 1–80 (2003)

<sup>12</sup> J. Oxley et al “Raman and infrared fingerprint spectroscopy of Peroxide-based explosives,” Applied Spectroscopy Vol 62:8, 906-915 (2008)